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

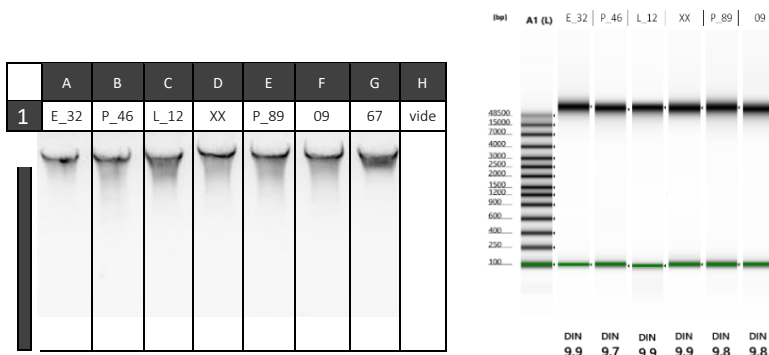
Never send the entire DNA stock to the PGTB, only aliquots.

Sample preparation

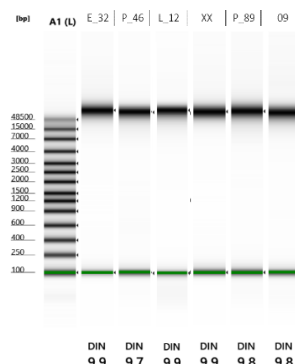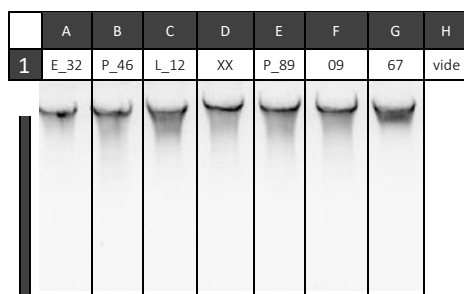- ⇒ Quantify DNA by **fluorescence**, then dilute it between 20 and 30 ng/μL. If concentrations are < 10 ng/μL, inform the PGTB. Once samples have been diluted, check the concentration again using fluorescence.
- ⇒ Check the integrity of DNA using an agarose gel or, ideally, a capillary electrophoresis device (Bioanalyzer, Tapestation, etc.). Annotate the gel, specifying the name of samples or their position on the plate.

From 1 to 16 samples:

- ⇒ Pipet (at least) **30 μL** of the diluted DNA into **LoBind Safe-Lock tubes**.
- ⇒ Identify tubes with **the acronym of the project declared in the project form** and the sample name. Close and seal the tubes with Parafilm and place them in a box also bearing the project acronym.

For more than 16 samples:

- ⇒ Use **96-well PCR plates that are transparent, skirted and non-break-away**. Identify each plate with **the acronym of the project declared in the project form** and a plate number: PROJECT X – PLATE 1.
- ⇒ Pipet (at least) **30 μL** of diluted DNA into the plates, **leaving one well empty without water or DNA on each plate as follows**: plate 1 = H1 empty, plate 2 = H2 empty, ..., plate 12 = H12 empty.
- ⇒ Heat seal the plates with suitable films (be careful not to deform them by overheating) or seal the plates with compatible cap strips. If you have used caps, wrap the plates in Parafilm to keep the caps firmly in place.

Depositing or sending samples

- ⇒ Complete [the sample submission form](#) to deposit samples at the PGTB or to send them by courier. Parcels are received from 9:00 a.m. to 5:30 p.m., Monday to Friday, at the following address:

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- ⇒ For shipping by courier, place ice packs at the bottom and sides of the package. Place the box containing the tubes or plates in the centre (for plates, separating them with cardboard to prevent them from being crushed). Fill any empty spaces with bubble wrap or cardboard so that the plates are securely held in place in the package.
- ⇒ PGTB will notify you by email when your samples have been received. **Any samples that do not meet the conditions specified in the email or in these requirements cannot be processed within the stated time frame. An additional charge may be applied or a new sample shipment may be requested.**

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⇒ Send the list of sample name to pgtb-lab@inrae.fr.

For more than 16 samples:

- ⇒ Once we have received sample submission form, you will receive an email from PGTB with a link to directly access the file for submitting plate designs for your project on our RESANA platform.
- ⇒ Click on the link and the plate plan submission file will open in your web browser. Complete the plate plans following the instructions provided: **only letters/numbers/hyphens are permitted**.
- ⇒ You can copy/paste the plate designs using the keyboard shortcut 'Ctrl + V'.
- ⇒ **The file is automatically saved, so you do not need to do anything**. Once all the plate designs have been completed, close the file by clicking on the cross in the top right-hand corner.

	Rédaction	Zoé Delporte - AI	
	Approbation	Erwan Guichoux - IR	