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|   |   |                             |
|---|---|-----------------------------|
|  | <b>Requirements for sending samples for<br/>Illumina Sequencing (ready-to-load)</b><br>(French version on page 1) | <b>Instruction</b>          |
|   |   | <b>Version : 28.04.2026</b> |
|   |   | <b>Pages : 2/2</b>          |

**Never send the entire DNA stock to the PGTB, only aliquots.**

### Pool preparation

- ⇒ After preparing your pool, check **its concentration, ideally** by qPCR, otherwise by fluorescence, as well as **its size**.
- ⇒ Place an aliquot of (at least) **30 µL** of the **5nM** pool in a 1.5mL or 2mL Eppendorf LoBind tube, close and seal the tube with Parafilm.
- ⇒ Identify the tube with a label indicating the project acronym and a unique number if there are several tubes.

### Sample information in the pool

- ⇒ Use only **unique** identifiers for your samples, controls and replicates, and use only letters, numbers and hyphens, with no spaces or dots. Identifiers must not exceed 50 characters.

- ⇒ Prepare a summary file containing the following information for each sample:

- Unique identifiers and associated indices (1 index = 1 sample)
- Number of cycles per read:
  - *Read1Cycles* : 50
  - *Index1Cycles* : 8
  - *Index2Cycles* : 8
  - *Read1Cycles* : 50
- Loading concentration : 750pM
- PhiX percentage : 5%

| ID    | Index i7   | Index i5   |
|-------|------------|------------|
| E32   | GACCAAGTTA | GATTCACGAC |
| P-46  | CTCCGCTTAT | CGTAGAACAG |
| L12   | ATTAAGTTCG | ACCTAAGAGC |
| XX    | GTGGCAACAT | ATGGCTTCTA |
| P-89  | TGACCTGCTG | TACCTAAGTG |
| 9B    | CGCAGCTAGA | CCTTACCTAT |
| 67    | ACGCTAATCT | ATATAAGGCG |
| Alpha | GTTGTTACCT | AGCATCGTGC |

**Please note:** In a **single** sequencing run, indexes must ALL be of the same size and format (dual/single).

**Note:** If you do not specify a loading concentration, we will load the cassette according to Illumina recommendations with a library concentration of 750pM and a PhiX concentration of 5%.

- ⇒ Send this file with sample <-> index matches to [pgtb-lab@inrae.fr](mailto:pgtb-lab@inrae.fr) by e-mail, specifying the **concentration** and **average pool size**, as well as the **references of the library and index kits** used to prepare the libraries.

### Depositing or sending samples

- ⇒ Complete [the sample submission form](#) to deposit samples at the PGTB or to send them by courier. Parcels are received from 9:00 a.m. to 5:30 p.m., Monday to Friday, at the following address:

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- ⇒ For shipping by courier, place ice packs at the bottom and sides of the package. Place the box containing the tubes. Fill any empty spaces with bubble wrap or cardboard so that the plates are securely held in place in the package.
- ⇒ PGTB will notify you by email when your samples have been received. **Any samples that do not meet the conditions specified in the email or in these requirements cannot be processed within the stated time frame. An additional charge may be applied or a new sample shipment may be requested.**

|   |             |                     |   |
|---|-------------|---------------------|---|
|  | Rédaction   | Zoé Delporte - AI   |  |
|   | Approbation | Erwan Guichoux - IR |   |